

## Firmware Update Notification

 Urgent Action Required   
 Information Only 

<b>NEW FIRMWARE VERSION</b>	<b>EFFECTIVE DATE: NOVEMBER 2023</b>										
<b>TC LAB REVISIONS A - C</b> TC5 LAB, TC10 LAB, TC15 LAB → Firmware Version 1.99	<b>TC LAB REVISION D</b> TC10 LAB, TC15 LAB → Firmware Version 2.01										
<b>UPDATED COMMAND - CONST:THERM</b>											
<p><b>Subject: TC LAB Firmware Update</b></p> <p>The TC LAB Instrument firmware has received an update to include the ability to manually select a fixed and persistent bias current for a thermistor-type sensor. This update is tied to the remote command <i>CONST:THERM</i> and allows for an auto-ranging bias current or a fixed and persistent bias current when creating a custom sensor. The selected bias current will be automatically recalled and set to the active bias current from memory by the firmware in the event that the TC LAB is power cycled from the front or rear panel or if the sensor load is removed/disconnected.</p> <p><b>CHANGES AND FEATURES INCLUDE:</b></p> <ul style="list-style-type: none"> <li>Fixed bias currents available for custom thermistor</li> </ul> <p><b>REMOTE COMMAND - CONST:THERM</b> (see page 63 of the LAB Series <a href="#">Command Set</a>)</p> <p>To create a sensor with a fixed bias current, append a suffix to the name of the sensor when it is created. The format of the suffix is:</p> <p style="padding-left: 40px;">&lt;sensor name&gt;.Fn</p> <p>where &lt;sensor name&gt; is the symbolic name of the new sensor, and n indicates the bias current value, as would be passed to the <i>TEC:BIAS</i> command. For the <i>TEC:BIAS</i> command, the inputs are shown in Table 1.</p> <table border="1" data-bbox="1263 961 1498 1199"> <caption>Table 1</caption> <thead> <tr> <th>INPUT</th> <th>BIAS CURRENT</th> </tr> </thead> <tbody> <tr> <td>1</td> <td>10 μA</td> </tr> <tr> <td>2</td> <td>100 μA</td> </tr> <tr> <td>3</td> <td>1 mA</td> </tr> <tr> <td>4</td> <td>10 mA</td> </tr> </tbody> </table> <p>To create a sensor that will used a fixed bias, use <i>CONST:THERM</i> remote command. For example, the command to create a 50kΩ @ 25°C thermistor with a fixed 100μA bias current is:</p> <p style="padding-left: 40px;">CONST:THERM My50k.F2, 9.6911E-04, 2.1014E-04, 8.8019E-08</p> <p>This creates a new sensor named “My50k” that, when selected, will set a fixed and persistent 100μA bias current. Once created, the custom sensor can be selected from the front panel touchscreen. Additionally, the <i>TEC:SENSOR</i> command can be used to select this sensor.</p> <p style="text-align: center;"><b>REVIEW THE DATASHEET AND COMMAND SET FOR DETAILED INFORMATION ON THE UPDATED CUSTOM SENSOR SPECIFICATIONS AND OPERATION.</b></p> <p><b>Action Required:</b> Update Firmware Version - follow updating firmware instructions on page 19 of the TC LAB Instrument Datasheet found here: <a href="http://www.teamwavelength.com/download/Datasheets/tc_lab.pdf">www.teamwavelength.com/download/Datasheets/tc_lab.pdf</a></p> <p style="text-align: center;"><b>NOTE: INSTALLING THE WRONG REVISION OF FIRMWARE WILL REQUIRE RECALIBRATION AT THE FACTORY.</b></p>		INPUT	BIAS CURRENT	1	10 μA	2	100 μA	3	1 mA	4	10 mA
INPUT	BIAS CURRENT										
1	10 μA										
2	100 μA										
3	1 mA										
4	10 mA										
Please contact Sales at 406-587-4910 or email <a href="mailto:sales@teamwavelength.com">sales@teamwavelength.com</a> for details											